


# Time Lapse

## Note

Use this time lapse mode when a time interval of a second or more between images is needed.

1. Open **QCapture-Pro** or **Image-Pro**
2. Select the **Acquire** file menu then select the **Video/Digital** option. In the camera control window you will find the **Image** tab displayed on the right (you may need to click on **More>>>** at the bottom of the window to reveal this)
3. Tick **Enable Multiple Image Capture**
4. Type or select the desired number of images (100 in our example)
5. Click **Set Time** and enter a time interval between images in the **Time Interval** dialog
6. Select the **File** option and click **Browse** to select a save location
7. Click **Snap** to grab a series of images.

**Note** If the **Set Time** button is greyed out, deselect the **Use minimum possible time interval** option.

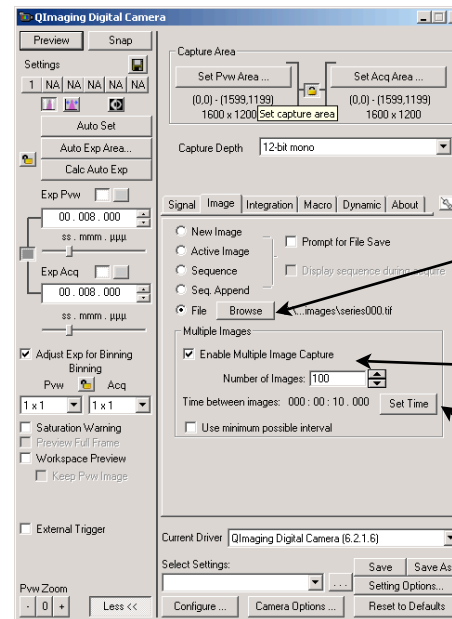

File

Enable Multiple Image Capture

Set Time

